4-Channel Active Base for the Hamamatsu S13361-3050AE-08 Rev. P1

Features

Supports the Hamamatsu S13361-3050AE-08 8x8 array of 3mm MPPCs

"HFN" Variant: Horizontal signal connector on the front, array located on the front, narrow format

3-side tileable installation

Four encoded position signals for event centroid calculations: X+, X-, Y+, Y-

DC-coupled signal path

Low power consumption

Patented diode-coupled charge division readout, superior to traditional resistive readout


Improved spatial uniformity


Faster rise time


Reduced image noise

Precision temperature sensor

Four mounting holes for #4 or M3 hardware

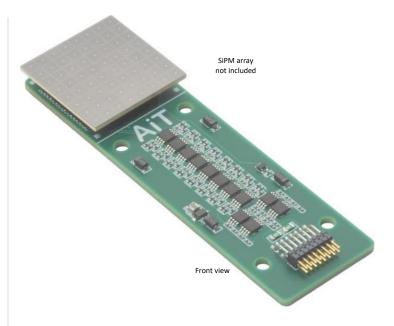

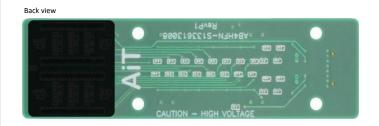

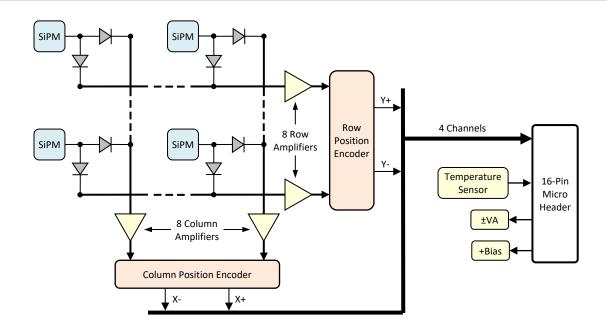

AB4HFN-S133613008

Specifications

Position Signal Outputs

Encoding Charge division multiplexed to

4 output channels: X+, X-, Y+, Y-

Gain 750Ω transimpedance gain

Output voltage $0 \rightarrow -1V$ into 100Ω load

Output impedance 100Ω

50mA maximum Output current

Temperature Sensor

Output voltage 500mV + 10mV per °C

Output current 10mA Output impedance 100Ω ±0.5°C Accuracy

Bias Voltage +56V typical (refer to SiPM data)

Voltage clamp 68V Zener diode

375mW maximum

Amplifier Voltage (±VA) ±2.8V → ±5.5V DC maximum

Current ±40mA typical

(Iq, no signal, no load)


Signal Connector Horizontal 16-pin 2-row header

0.050" pin pitch


Samtec FFSD-08-D-XX.XX-01-N Mating assembly

(XX.XX = length in inches)

Channel Map

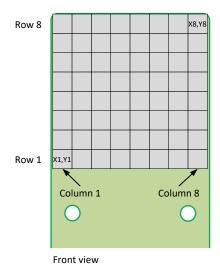

Signal Connector

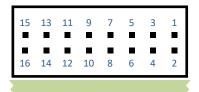

PCB Side View

| Pin | Function | Pin | Function |
|-----|-------------|-----|----------|
| 1 | Temperature | 2 | Ground |
| 3 | X- | 4 | Ground |
| 5 | X+ | 6 | Ground |
| 7 | -VA | 8 | Ground |
| 9 | +VA | 10 | Ground |
| 11 | Υ- | 12 | Ground |
| 13 | Y+ | 14 | Ground |
| 15 | +Bias | 16 | Ground |

4-Channel Position Encoder

Row and Column Encoder Weights

| Row# or Col# (for X- or Y-) | Row# or Col# (for X+ or Y+) | Fraction (ideal) | Fraction (actual) | % Error | Notes |
|--------------------------------|--------------------------------|---------------------|----------------------|---------|--------------------------------|
| 1 | 8 | 0.1250 | 0.1250 | 0.00 % | |
| 2 | 7 | 0.1230 | 0.1230 | -0.68 % | • |
| 3 | 6 | 0.2300 | 0.2463 | -0.51 % | Sum of X- and X+ fractions |
| 4 | 5 | 0.5000 | 0.5000 | 0.00 % | or Y- and Y+ fractions |
| 5 | 4 | 0.6250 | 0.6250 | 0.00 % | = 1.1250 |
| 6 | 3 | 0.0230 | 0.0230 | 0.00 % | Independent of signal position |
| 7 | 2 | 0.7300 | 0.7300 | 1.44 % | macpendent of signal position |
| 8 | 1 | | | | |
| 8 | 1 | 1.0000 | 1.0000 | 0.00 % | |

Note: Errors exclude component tolerances

Output Signals

X- = (SiPM signal) * (encoder gain) * (X- fraction)

X+ = (SiPM signal) * (encoder gain) * (X+ fraction)

Y- = (SiPM signal) * (encoder gain) * (Y- fraction)

Y+ = (SiPM signal) * (encoder gain) * (Y+ fraction)

Typical event position calculation:

X column = (X + - X -) / (X + + X -)

Y row = (Y+ - Y-) / (Y+ + Y-)

Example

SiPM signal at column 4, row 3 (excluding encoder gain)


X- = (Column 4 signal) * 0.5000


X+ = (Column 4 signal) * 0.6250


Y- = (Row 3 signal) * 0.3731


Y+ = (Row 3 signal) * 0.7500

Typical Signals

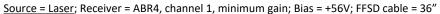


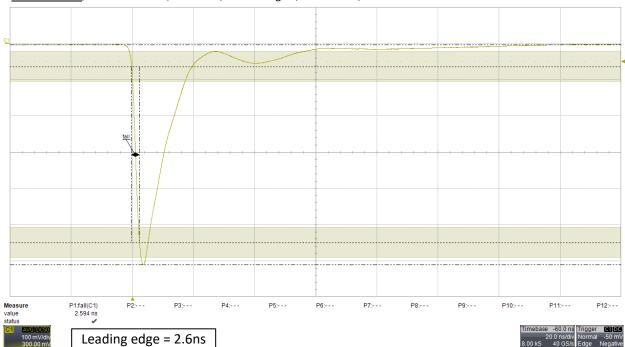

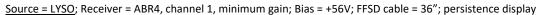


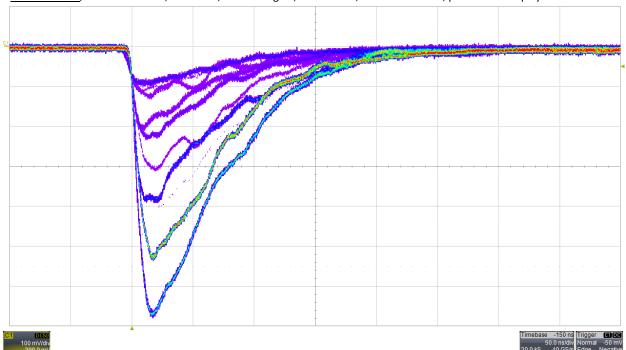


Mechanical

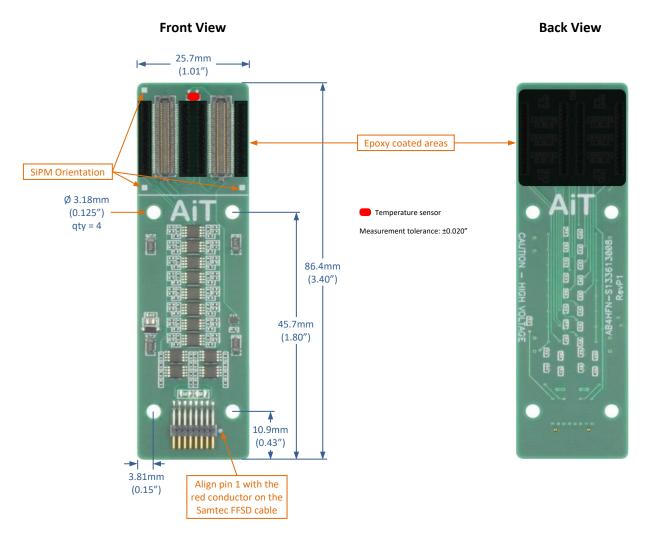

Array Location

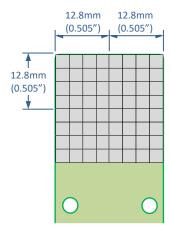

4-Channel Active Base Readout Kit

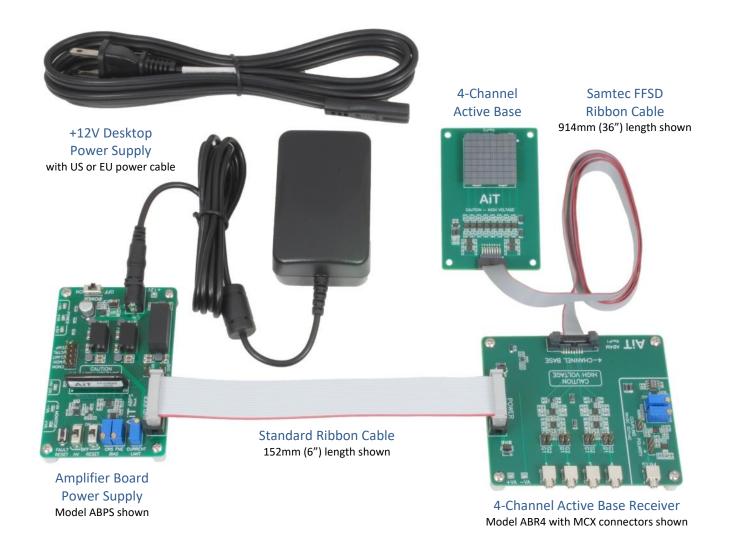

Components

Each component is available separately. Refer to each datasheet for details.

The Active Base includes a 914mm (36") Samtec FFSD micro-pitch ribbon cable.

The Amplifier Board Power Supply includes a 12V desktop power supply and a HV80 bias voltage power supply.

The 4-channel Active Base Receiver includes a 152mm (6") power supply ribbon cable and a breakout board to connect any external power supply.

AB4HFN-S133613008

Safety Information



WARNING – High Voltage

- High voltage may be present during operation
- High voltage stored on capacitors may be present after power is removed
- Improper handling may result in personnel injury or equipment damage

This high-voltage device must be used only by personnel trained and qualified in safe handling, installation, and operation of high-voltage equipment.



CAUTION – Electrostatic Discharge (ESD) Sensitivity

The circuit board can be damaged by electrostatic discharge. Observe precautions for handling electrostatic sensitive devices. Handle only at static-safe workstations.

High-Gain Photodetectors

High-gain photodetectors such as silicon photomultipliers may conduct damaging currents if exposed to high optical signal levels while the bias voltage is applied, or if the bias voltage exceeds the recommended operating range. These devices must be operated only in low-light conditions, and only within the manufacturer's recommended bias voltage range.

Handling and Disassembly

This product may be provided with a protective enclosure. Disassembled enclosure components and circuit boards may contain sharp edges. Take appropriate safety precautions while assembling or disassembling the enclosure and handling disassembled components.

Indoor Use Only

Do not operate this product in a wet or damp environment. Do not operate in an explosive atmosphere.

Use of this product, and AiT Instruments' liability related to use of this product, is further governed by AiT Instruments' standard terms and conditions of sale, which were provided upon purchase of this product.